



## **BUNNY SNUGGLE**

Congratulations on the purchase of your new BUNNY SNUGGLE the safest form of rabbit restraint. Please assure you have purchased the correct size for your rabbits. You are now ready to start your first stress free rabbit procedure.

First open your SNUGGLE on a flat surface with the rear extension facing away, see #1.

With your subjects head tucked under your arm place the subject on the SNUGGLE facing #8.

While holding the subject firmly raise panel #1 and hold with the thumb and forefinger of your hand. Close flap #2 over #1 then proceed to cover #2 with #3, assure a firm closure. Now fasten #5 over #4 behind the neck. The eyes should now be covered with #7 secured over #6. This will keep the subject calm during procedures.

Panels #7 and #4 as well as #6 and #5 may be interfaced for a more secure hold of the head (gavage, eye movement etc.)

The SNUGGLE may be used in many configurations for various testing and sampling procedures. Should you have any suggestions or need more information please contact us.

UNO Roestvaststaal BV  
P.O. Box 15  
6900 AA ZEVENAAR  
The Netherlands  
Phone: +31 316 524451  
Fax.: +31 316523785  
Email: [info@unobv.com](mailto:info@unobv.com)